


## Xdrop® Sort Idle function

- The new Xdrop Sort Idle function that is not described in Xdrop Sort Manual v 4.0.
- This function allows you to pause sorting on an individual lane. Using it has no impact on other lanes. You can use it on multiple lanes.
- You must activate the Idle function if no sorting events occur on a given lane within 20 min of the start of the run to prevent instrument damage.
- **Idle** pauses sorting on one lane only. **Pause Sorting** pauses sorting on all active lanes.

### How to use the Xdrop Sort Idle function

1. Using the buttons along the bottom of the Xdrop Sort screen to select a lane. Lane 1 is selected in the screenshot below.
2. The **Idle** button is at the right of the screen at the top of the button stack. Press it to activate the Idle function. Sorting on the displayed lane will pause.
3. Press the button again to restart sorting on that lane.

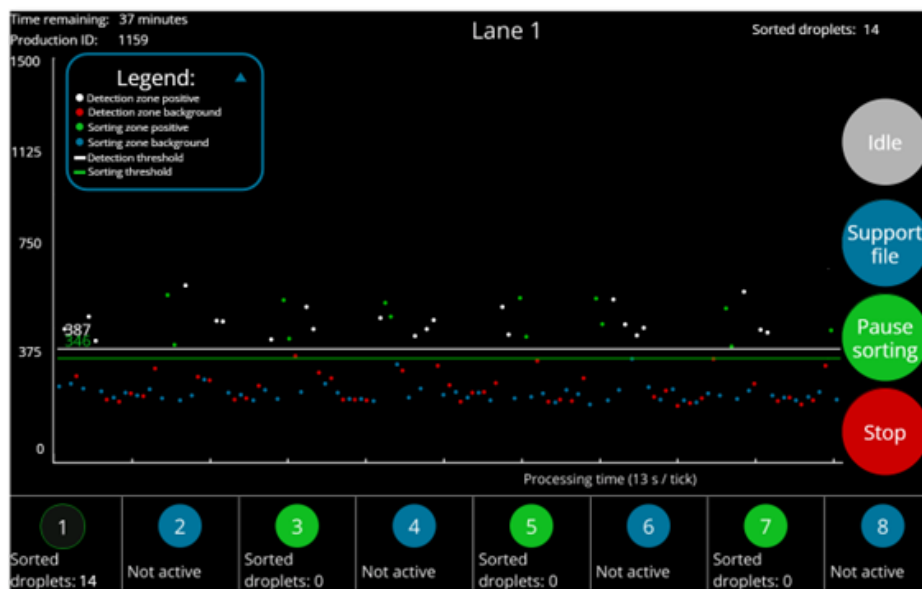

Please refer to the Xdrop Sort Manual version 4.0 or later for more information.