

## How to update xdrop™ Instrument's Software

The software version of the Xdrop™ instrument is reported on the bottom-left corner of the instrument screen.









### Software Update

1. Download the updated software from our support website [here](#). Go to the end of the page and download the Software update zip file:

<p><b>Xdrop™ Software Downloads</b></p> <p>Here you will find the latest versions of Xdrop™ Software for downloading and updating your Xdrop™ instrument.</p>	<p><b>Latest Updates</b></p> <p>Xdrop™ Software Update. Version 1.1.0. Released 26th November 2019</p> <p><a href="#">Download version 1.1.0</a></p>
---	--

Note that most recent version might be different from the one reported in the above screenshot.

2. Unzip the downloaded file and save the unpacked files on the USB dongle provided with the instrument. This is the folder structure at the root of the USB dongle, after files have been successfully unzipped to the root of the device:

 colibri_imx6	File folder	
 flash_blk	Disc Image File	1 KB
 flash_eth	Disc Image File	1 KB
 flash_mmc	Disc Image File	1 KB
 software_version	Text Document	0 KB
 xdrop_update	Text Document	0 KB

3. Plug the USB dongle into the back of the instrument. A new button stating “Software Update” should appear on the instrument screen.



4. Press “Software Update”. Wait for the update to finish and then remove the dongle. **Note:** Screen will go black during the first part of the update and will eventually re-boot to complete the update.

If you need guidance throughout the process, get in touch with us at [support@samplix.com](mailto:support@samplix.com).